# LDZM 系列

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Vertical Pressure Steam  
Sterilizer

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型号	LDZM-40L	LDZM-60L	LDZM-80L
	LDZM-40L-I	LDZM-60L-I	LDZM-80L-I
容积	40 立升	60 立升	80 立升

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### I. Basic structure

1. The shell case is made of high temperature thermoplastic material with outstanding impact resistance and stainless steel.
2. The pot body of sterilizer is made of food grade stainless steel (304).
3. The lid opening and closing device adopts the lever type multi-link synchronous telescopic structure, which makes the lid and the pot body open and close flexible and light, safe and reliable.

### II、 Control and operation

The console uses the LCD window graphic dynamic display, the function settings are all selected by scrolling graphic and text, the whole sterilization phase is monitored by micro-computer graphic display and automatic cycle control, and the automatic shutdown is initiated when sterilization phase ends (with an alarm).

### III、 Function description

1. It has a one-button sterilization function selection, which can make a quick and clear choice based on different items.
2. With air-cooled chilling device to ensure that the pot body is quickly cooled at the end of sterilization, thus reducing the time to open the lid.
3. It has the function of preheating and melting of sterilized items and has time and temperature selection for the items to be dissolved.
4. A USB interface is available for the convenience of monitoring process with connection to a computer and of collecting data.
5. Sterilization reservation function is available so that users can start it at a set time.
6. Automatic fault detection and judgment system is in place.
7. The full-automatic control model has the function of discharging the low temperature air and exhausting steam when sterilization procedure ends.
8. Add print function, print the current temperature and sterilization time.

### IV、 Safety devices

1. It has a safety interlocking device, structure of which is featured by electronic and mechanical interaction, thus ensuring that it automatically locks the lid when there is pressure in the pot and avoid any danger resulted from improper operations.
2. Double protection system with the mechanical safety relief valve and electronically controlled overvoltage protective device.
3. It has heater anti-drying protection function and leakage protection system.

## Technical Parameter

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参数 / 型号 Parameter / Model	LDZM-40L LDZM-60L LDZM-80L	LDZM-40L-I LDZM-60L-I LDZM-80L-I
配置 Configuration	自控型 Semi-automatic control	全自控型 Full-automatic control
主体材质 The main material	不锈钢 Stainless steel	不锈钢 Stainless steel
容积 / 毛重 Volume / GW	93/95.5/110Kg(40L/60L/80L)	97/99/113.5Kg(40L/60L/80L)
电压 / 功率 Voltage / Power	220V / 3.5Kw	220V / 3.5 Kw
包装尺寸 (长 x 宽 x 高 mm) Packing size (L X W X H mm)	780×660×1160 (40L/60L)	780×660×1320 (80L)
产品尺寸 (长 x 宽 x 高 mm) Product size (L X W X H mm)	570 X 690 X 1020 (40L) 570 X 690 X 1020 (60L)	570 X 690 X 1170 (80L)
容器容积 (直径 x 高度 mm) Sterilization chamber volume(D X H mm)	370 X 380 (40L) 370 X 560 (60L)	370 X 750 (80L)
网篮尺寸 (直径 x 高度 mm) Basket Size(D X H mm)	350 X 190 (40L) 350 X 360 (60L)	350 X 190 & 350 X 360 (80L)
时间设定范围 Time setting range	0-99 h	0-99 h
额定工作压力 Rated working pressure	0.217MPa	0.217MPa
灭菌温度设定范围 Sterilization temperature setting range	50~134°C	50~134°C